

IDEXX[®] In-House (IVLS)

User Guide



877-999-2383 | covetrus.com

© 2025 Covetrus, Inc. All rights reserved. No copying without consent.



TABLE OF CONTENTS

Overview	. 3
Advanced Options	3
Mapping	. 4
Select the Device	. 5
Place the Order	. 5
Notifications	. 6

covetrus 📚

Overview

The new IDEXX IVLS integration has the ability for Avimark to communicate with multiple IVLS devices in the practice. This document will go through the process of configuring the settings.

Advanced Options

Enter the IDEXX In-House credentials under Advanced Options in Avimark.

- 1. Click the gear icon () on the speedbar.
- 2. Search for 'idexx in-house'.
- 3. Set Active to True.
- 4. Enter the API Password and API Username.

5. Enter the LIMS ID.

The LIMS ID with username and password will let the program know what IVLS machines are in-house.

6. Click Done.

Coptions Maintenance			\times
<u>File E</u> dit <u>V</u> iew <u>H</u> elp			
Search idexx in-house			
Options Lab Instruments DEXX In-House Active API Password API Username LIMS ID			
Values Security			
Default Value	Change		
		 Do	ne

covetrus 📚

Mapping

Because the IDEXX IVLS In-House is a new integration, species and breed mapping will be performed automatically. However, mapping can also be done manually, if needed.

- 1. From the menu, click on Work with > System Tables.
- 2. Select the **Species** or **Breed** table.
- 3. Double-click on the species or breed that requires mapping.
- 4. Click on the Mapping tab.
- 5. Under **Code** next to IDEXX In-House, double-click to get the list and select the mapping code for the species or breed selected.
- 6. Click OK.

😤 Change Breed Entry - Akita Mix		×
Table Breed Table Details Mapping		
Integration	Code	
IDEXX Reference Laboratories	AKITA	
IDEXX In-House	N	
	6	
*** - means a mapping proposal		
0	DK	Cancel



Select the Device

When there are multiple IVLS machines, a prompt will appear for the user to select the desired device for sending the lab request. Avimark will remember the last IVLS selected on the computer but a different device can be selected from the drop-down list.

Note: This prompt will not appear if there is only one IVLS.

- 1. In the Patient area, right-click > Create Requisition.
- 2. Select the IVLS device from the drop-down list.

Please select device		×
Please select the appropria	te /VLS device	
IVLS	PTH999900000627 - IVLS	-
	PTH999900000625 - IVLS	
	PTH999900000626 -LALS	
	PTH999900000627 - IVLS	
	PTH999900000628 - IVLS	

3. Proceed with placing the order.

Place the Order

After the IVLS device is selected, the screen will appear for the user to click the button to place the order.

integration.vetconnectplus.com/ui?token=5efa0cdf01ff7c6be0ca9077e0640d0a&corderId=100264964	-		×
IDEXX DIAGNOSTICS			
FLUFFY L'HEUREUX			
IDEXX VetLab Station Pre-Anesthetic Panel (PreAn)			
* Veterinarian			
Nikki L'Heureux, DVM			
Search All Tests		Ord	er



Notifications

IDEXX In-House lab results, when complete, will appear under Notifications.

- **1.** Right-click on a completed result.
- 2. Click on Import Results.

9		Notifications	
Time 🔻	Category	Message	
9:18 AM	IDEXX In-H	L'Heureux Eluffy - results	ready (F)
9:18 AM	IDEXX In-H	L'Het Acknowledge	ady (F)
9:08 AM	Zoetis Con	L'Het Jenned Provide	TED
9:08 AM	Zoetis Con	Dane	
9:08 AM	Zoetis Con	Claus, Snowy - COMPLE	TED
9:08 AM	Zoetis Con	Conner, Day Day - COMF	PLETED
9:08 AM	Zoetis Con	Bradley, Shaggy - COMP	LETED
9:08 AM	Zoetis Con	Carswell, Muffin - COMPL	ETED
9:08 AM	Zoetis Con	Danes, Fluffy - COMPLE	TED
9:08 AM	Zoetis Con	Danes, Jerry - COMPLET	ED

3. Requisition Management will load the IDEXX In-House labs and download the results.

File Edit View Help								
		<u>File Edit View</u>	Help					
Requisitions Download Requisitions		Requisitions I	lownload Requisitions			_		
Lab Type IDEXX In-House - Loading		Lab Type IDE	EXX In-House 👻	Load				1
Created Date 🤍 Lab Client Name Lab Patient Name Result Status RequisitionID Patient	t AccessionI	Created Date 04-30-25 9:17a 04-30-25 9:15a	♥ Lab Client Name	Lab Patient Name	Result Status Final Final	RequisitionID 100264965 100264964	Patient Fluffy Fluffy	Accessio
		04-22-25 8:30a 03-25-25 2:18p	Tim Grodon Morehead	Felix Keika	Final Final	100263979 100258631		
<no data="" display="" to=""></no>								